

CesiumTaq LA

Cat #: 210



Amount: 100 µl (2000 x 25 µl reactions)

Shipping conditions: Ambient temperature

Storage conditions: -20°C

Thermostability: Retains at least 85% activity after 1 hour at 95°C

Expiration: On tube label

PRODUCT DESCRIPTION:

CesiumTaq LA is a DNA polymerase mixture containing CesiumTaq, a double cold-sensitive mutant of Taq DNA polymerase. Due to its suppressed activity at low temperatures this enzyme is designed for hot-start PCR performance. The Long-Accurate feature allows for amplification of longer products with higher fidelity and accuracy. LA enzymes are not recommended for use with dUTP. 10x buffer composition is: 500 mM Tris-Cl pH 8.3, 160 mM ammonium sulfate, 0.25% Brij 58, and 25 mM magnesium chloride.

TYPICAL PCR PROTOCOL for a 25 µl reaction:

Reagent	Volume	Final Concentration
10x Taq Mutant Reaction Buffer	2.5 µl	1x
dNTP mix (10 mM each)	0.5 µl	200 µM each
Left Primer	variable	200 nM
Right Primer	variable	200 nM
DNA template†	variable	0.1-100 ng
PCR Enhancer Cocktail (recommended)*	12.5 µl	1X
CesiumTaq LA	0.05-.25 µl**	
De-ionized distilled H2O	Adjust final volume to 25 µl	

† DNA amount depends mostly on genome size and target gene copy number.

* For optimal performance, we recommend using one of our PCR Enhancer Cocktails (PEC-1, PEC-1GC, PEC-2, or PEC-2-GC) or 1.3M Betaine, a general PCR enhancer.

** To determine specific optimal enzyme concentration, we strongly recommend an enzyme titration test for each target. A good starting amount of the enzyme per 25 µl reaction is 0.05 µl. Targets larger than 1 kb may require more enzyme.

CYCLING CONDITIONS:

1. Denaturing: 94° for 2-8 minutes for 1 cycle
2. Denaturing: 94° for 40-60 seconds
3. Annealing: 50°-68° depending on the specific T_m primers for 40-60 seconds
4. Extension: 68° for 2 min/kb target
5. Repeat steps 2-4 for 25-40 cycles

Please visit us on the web at www.klentag.com for troubleshooting and detailed protocols.

REFERENCES:

Kermekchiev, M.B., et al. (2003) Cold-sensitive mutants of Taq DNA polymerase provide a hot start for PCR. Nucl Acids Res. 31, 6139-6147.