

## 5x Klentaq LA PCR Kit Cat #: 130

**Amount:** 5 X 0.5 ml (500 x 25 ul reactions)

**Shipping conditions:** Ice pack

**Storage conditions:** for best performance, store at -20°C

**Shelf life:** At least 1 year if stored at -20°C and 10 freeze/thaws or at least 3 months if stored at 4°C.

### PRODUCT DESCRIPTION:

Our 5x ready-to-use PCR kit contains Klentaq LA, the Long Accurate version of Klentaq1 DNA Polymerase, a 5'-exonuclease deficient Taq polymerase (an N-terminal deletion of Taq) with improved fidelity and thermostability. This kit can be used for regular, as well as real-time PCR. It contains everything necessary for a PCR reaction to work perfectly, just add your template, primers/probes and water. For real-time reactions you may need to add a fluorescent dye as an alternative to probes. 5X composition is: 5x KlentaqLA DNA Polymerase, 1 mM dNTPs, 250 mM Tris-Cl pH 9.2, 80 mM ammonium sulfate, 0.125% Brij 58, and 17.5 mM magnesium chloride.

### TYPICAL PCR PROTOCOL for a 25 µl reaction:

Reagent	Volume	Final Concentration
5x KlentaqLA PCR Kit reagent	5 µl	1x
Left Primer	variable	0.2 µM
Right Primer	variable	0.2 µM
DNA template <sup>†</sup>	Variable	0.5-100ng
Betaine 5M*	6.5µl (optional)	1.3 M
de-ionized distilled H <sub>2</sub> O	Adjust final volume to 25 µl	-

<sup>†</sup> DNA amount depends mostly on genome size and target gene copy number.

\*Betaine is a general PCR enhancer. It usually improves the yield and specificity of amplification especially for longer targets.

### CYCLING CONDITIONS:

1. Pre-incubation: 94° for 2 minutes for 1 cycle
2. Denaturing: 94° for 40-60 seconds
3. Annealing: 55°-70° depending on the specific primers (5° less than T<sub>m</sub>) for 40-60 seconds
4. Extension: 68° for 2 min / 1kb target
5. Repeat steps 2-4 for 25-40 cycles

### REFERENCES:

Barnes, W.M. (1994) PCR amplification of up to 35 kb DNA with high fidelity and high yield from bacteriophage templates, PNAS 91, 2216-2220.

**Please visit us on the web at [www.klentaq.com](http://www.klentaq.com) for troubleshooting and detailed protocols.**

### Notice to Purchaser

DNA Polymerase Technology products may not be resold, modified for resale or used to manufacture products without an agreement with DNA Polymerase Technology, Inc. Klentaq1 by DNA Polymerase Technology is licensed under US Patent No. 5,436,149. No license for Klentaq LA to be used in a Polymerase Chain Reaction has been purchased by DNA Polymerase Technology, Inc.