

5x Cesium Taq PCR Kit

Cat #: 220



Amount: 125 µl enzyme and two 1.25 ml tubes of 5x Mastermix
(sufficient for 500 x 25 µl reactions)

Shipping conditions: Ice Pack

Storage conditions: -20°C F

Thermostability: Retains at least 85% activity after 1 hour at 95°C

Shelf life: At least 1 year from date of receipt under proper storage conditions.

PRODUCT DESCRIPTION:

Our 5x ready-to-use PCR kit contains CesiumTaq, a cold-sensitive double mutant of Taq polymerase. Due to its suppressed activity at low temperatures this enzyme is designed for hot-start PCR performance. This kit can be used for conventional, as well as real-time PCR. For real time applications you may need to add a fluorescent dye as an alternative to probes.

5x Mastermix composition is: 250 mM Tris-Cl, 80 mM ammonium sulfate, 0.13% Brij 58, 12.5 mM Magnesium Chloride, and 1 mM each dNTP. Final pH is 9.1.

TYPICAL PCR PROTOCOL for a 25 ul reaction:

Reagent	Volume	Final Concentration
5x Mastermix	5 µl	1x
Left Primer	variable	200 nM
Right Primer	variable	200 nM
DNA template†	variable	0.1-100 ng
PCR Enhancer Cocktail (recommended)*	12.5 µl	1x
Cesium Taq	0.05 – 0.25 µl **	
De-ionized distilled H2O	Adjust final volume to 25 ul	

† DNA amount depends mostly on genome size and target gene copy number.

* For optimal performance, we recommend using one of our PCR Enhancer Cocktails (PEC-1, PEC-1GC, PEC-2, or PEC-2-GC) or 1.3 M Betaine which is a general PCR enhancer.

** To determine specific optimal enzyme concentration, we strongly recommend an enzyme titration test for each target. Targets larger than 1 kb may require more enzyme or may benefit from the LA version of the enzyme.

CYCLING CONDITIONS:

1. Denaturing: 94° for 2 minutes for 1 cycle *
2. Denaturing: 94° for 40-60 seconds
3. Annealing: 50°-68° depending on the specific Tm primers for 40-60 seconds
4. Extension: 68° for 2 min/kb target
5. Repeat steps 2-4 for 25-40 cycles

Please visit us on the web at www.klentaq.com for troubleshooting and detailed protocols.

REFERENCES:

Kermekchiev, M.B., et al. (2003) Cold-sensitive mutants of Taq DNA polymerase provide a hot start for PCR. Nucl Acids Res. 31, 6139-6147.

Kermekchiev, M.B. et al. (2009) Mutants of Taq DNA polymerase resistant to PCR inhibitors allow DNA amplification from whole blood and crude soil samples. Nucl. Acids Res., 37 (5):e40 E pub.

Notice to Purchaser: DNA Polymerase Technology products may not be resold, modified for resale or used to manufacture products without an agreement with DNA Polymerase Technology, Inc. The Omni mutant DNA Polymerases are trademarked and patented (US 7,462,475, and US patent pending). No license for OmniTaq 3, to be used in a Polymerase Chain Reaction, has been purchased by DNA Polymerase Technology, Inc.